Brazil’s Ministry of Agriculture, Cattle Raising And Supply (“Ministério da Agricultura, Pecuária e Abastecimento – MAPA”)

BUREAU OF AGRICULTURE AND CATTLE RAISING SAFETY (“Secretaria de Defesa Animal – SDA”)

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The Secretary of the Bureau of Agriculture and Cattle Raising Safety (“SDA”), exercising the prerogative conferred by subparagraph II of Section 15, of Annex I, of Decree 4.629 of March 21, 2003, considering Directive Number 62 of August 26, 2003, and the content of Process number 21000.002596/2004-94, decides to:

Section 1 – Oficialize the validation of the methodology utilized in the pathogen detection system – A-BAX® for the detection of Salmonella spp in food samples, water and environmental samples – swabs as an alternative method equivalent to the reference method of the Ministry of Agriculture, Cattle Raising And Supply, according to the Anex of this Directive.

Section 2 – Sub-delegate the authority to the Director of the Department of Animal Safety (“Departamento de Defesa Animal”), to enact complementary directives necessary to the validation of reference methodologies adopted by the Ministry, in accordance to proposals presented by the Animal Laboratory Coordination (“Coordenação de Laboratório Animal”)

Sole Paragraph. The methodology addressed by this article shall be updated, whenever advisable by technological advances, through act of the Director of the Department of Animal Safety (“Departamento de Defesa Animal”), within the scope of his authority.

Section 3 – This Directive (“Instrução Normativa”) shall be in effect as of the date of its publication.

MAÇAO TADANO

ANNEX

VALIDATION OF THE METHODOLOGY UTILIZED IN THE PATHOGEN DETECTION SYSTEM – A-BAX® FOR THE DETECTION OF SALMONELLA SPP IN FOOD SAMPLES, WATER AND ENVIRONMENTAL SAMPLES – SWABS AS AN ALTERNATIVE METHOD EQUIVALENT TO THE REFERENCE METHOD OF THE MINISTRY OF AGRICULTURE, CATTLE RAISING AND SUPPLY (MINISTÉRIO DA AGRICULTURA, PECUÁRIA E ABASTECIMENTO – MAPA).

The work conducted had the objective of validating the methodology utilized by the A-BAX® System for the detection of Salmonella spp, in comparison with the methodology of reference utilized by the Ministry (Manual of Official Analytical Methods for Microbiologic Analysis for the Control of Animal originated products and Water – Directive number 62, of August 26, 2003) involving food and water samples and environmental swabs.

The work was initiated with an evaluation of the performance of the A-BAX® System benchmarked against the methodology of reference (Ministry), at LANAL – Microbiologia, through the utilization of homogeneous samples of 25g, prepared with meat matrix, fortified with a low concentration of the target microorganism (at a level close to that of the detection limit of the method of reference) and a high concentration of interferents. It was concluded that the A-BAX® System is comparable to the method of reference due to a statistical equivalence represented by a confidence level of 95%, based on the Chi square (x2) results, a level of correspondence =95%, sensitivity =98%, specificity =95%, false-negative occurrence =2.5%; false-positive occurrence =5%, for presenting a recovery of fractions equal to 92% in comparison with the method of reference.

70 different types of foods were simultaneously analyzed in four laboratories – LAPA/PE, LARA/SP, FOOD INTELLIGENCE/SP, and SFDK/SP, totalizing 1879 samples from the laboratories' routines. Out of the analyzed samples, 52 presented positive results and 1798 negative results in both methods; 28 samples presented positive results in the A-BAX® System and negative results in the method of reference, and one sample presented a positive result in the method of reference and a negative result in the A-BAX® System. These figures indicate that the A-BAX® System presents a sensitivity of 98.1%, a specificity of 98.5%, a false-negative level equal to 1.9%, a false-positive level equal to 1.5% and a relative precision equal to 98.5%.

Through the results obtains, it was concluded that the A-BAX® System is equivalent to the method of reference for the detection of Salmonella spp based on the results of Chi square (χ^2).

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